



Introducing the  
**ProFill *Vet***  
Veterinary  
Capsule Filler

*The ideal machine for: new products, small runs, & research.*

- Quickly fill hundreds of capsules
- Simple to operate
- Typical fill weights within 7-10% of mean
- Includes everything you need to begin filling capsules
- Easy cleanup
- Available in sizes #13 through #Su07

Item Number	Size Range	Number of Capsules per cycle	Output: capsules/hour
XV-VPFS A	Su7, 7, 10	25	250
XV-VPFS B	11	25	250
XV-VPFS C	12, 12el 13	42	420

Size	Outer Diameter (mm)	Locked Length (mm)	Volume	
			(oz)	(ml)
<b>Su07</b>	23.4	88.5	1 ¾	28
<b>7</b>	23.4	78.0	1 ½	24
<b>10</b>	23.4	64.0	1	18
<b>11</b>	20.9	47.5	½	10
<b>12el</b>	15.2	57.0	¾	7.5
<b>12</b>	15.3	40.5	¼	5
<b>13</b>	15.3	30.0	⅛	3

Available for **US\$ 595**

MANUFACTURED BY


**TORPAC**<sup>®</sup>

www.torpac.com info@torpac.com  
T: 1-973-244-1125 F: 1-973-244-1365  
333 Route 46, Fairfield, NJ 07004, USA

For more information go to:  
**www.torpac.com**

**Before you begin:**

Because each machine comes equipped to fill multiple size capsules, you must first install the proper Dome Plate Feet.

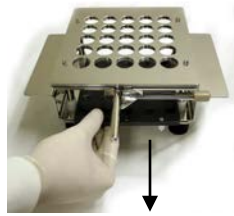

1. Remove the Dome Plate from the ProFill Vet as shown above.



2. Turn the Plate over and install the Dome Plate Feet that correspond with the capsule size to be filled.



3. Re-install the Dome Plate as shown above.



4. Load the capsules into the Caps Tray (longer narrower side first). Close the Locking Plate & Clamp to secure.



5. Before separating the capsules you must lock the bodies by, rotating the Cam Handle to desired degree (approx 60 to 80).



6. Separate the cap & body by pressing thumbs down on Caps Tray Handles and lifting the Caps Tray up with fingers.




7. Rotate the Cam Handle back to home position; the capsule bodies will fall down. Now the capsules are ready to fill.



8. Place the Powder Tray on the Filler by fitting the holes onto the Locating Pins on the Filler, keeping the Powder Pocket towards operator side. Pour pre-measured powder into filler.



9. Spread the powder evenly with the help of Powder Spreader. Push excess powder onto reservoir of Powder Tray.

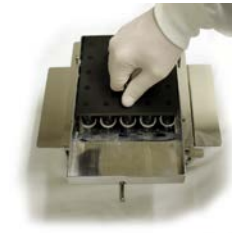

10. Tamp the powder of filled capsules with the help of Tamper as shown. Spread excess powder evenly and tamp again. Repeat this process until all capsules are completely filled.



11. After capsules are filled remove Powder Tray and place the Caps Tray back into position. **Prelock** the capsules by lifting the Lifting Plate as shown. **Do not try to lock capsules on Filler.**



12. Select corresponding Lock Spacer with capsule size being filled. Place selected Lock Spacer in the Locking Base as shown. This prevents over-locking of capsules.



13. Place the Capstray containing **prelocked** capsules on Locking Base by placing it over the Locating Pins.



14. Now lock the capsules by applying pressure with both hands (palms) as shown.



15. Remove capsules from the Caps Tray. You now have completely locked and filled capsules. With practice you can fill 400+ capsules per hour.

**TORPAC**<sup>®</sup>

www.torpac.com info@torpac.com  
T: 1-973-244-1125 F: 1-973-244-1365  
333 Route 46, Fairfield, NJ 07004, USA