







The Quickest Way to do Capsule Over-Encapsulation.


-  Simple to use.
-  Works with the ProFill 100 & ProFill OE Capsule Filling Systems.
-  Made of Acrylic and Stainless Steel.
-  Made for use with most capsule sizes.

The ProFill Capsule Filling System not included.


Capsule-in-Capsule Over-Encapsulation: To over-encapsulate a smaller size capsule into a larger capsule, you can use the ProFill 100 Orienter for the smaller size capsule, as it will orient the smaller size capsule even though the capsule is in the locked position. This attachment is available for immediate delivery. For size AAA only, this is a custom attachment, available in 6-8 weeks.

Capsule Size	Fits in DBcaps® size	Fits in Standard Capsule size
000		
00el		
00		000
0el		000, 00el
0	AAA, AAel	000, 00, 00el
1el	AAA, AAel	000, 00, 00el, 0
1	AAA, AAel	000, 00, 00el, 0, 0el
2el	AAA, AAel	000, 00, 00el, 0, 0el
2	AAA, AAel	000, 00, 00el, 0, 0el
3	AAA, AAel, AA, A	000, 00, 00el, 0, 0el, 1
4	AAA, AAel, AA, A	000, 00, 00el, 0, 0el, 1, 2
5	AAA, AAel, AA, A, B, D	000, 00, 00el, 0, 0el, 1, 2, 3


Using the Capsule-in-Capsule Over-Encapsulation Attachment.

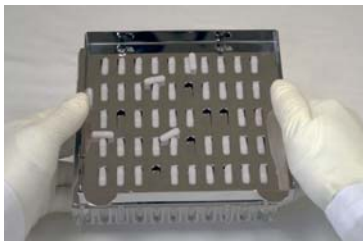

1. Pour 50+ trial capsules onto the Orienter. Shake back and forth moving them into the slots.

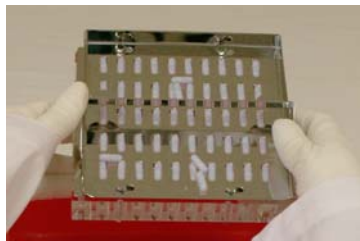

2. Lift gate and pour off excess trial capsules.

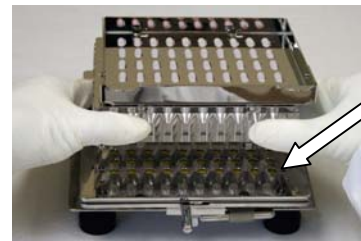


3. Place the Orienter locating feet into the holes further away from the operator on sheet "c".



4. Push the sliding portion of the Orienter to the left to drop trial capsules into the bodies of the outer capsule.



5. Remove the Orienter. Capsules are in every other row of the Filler.

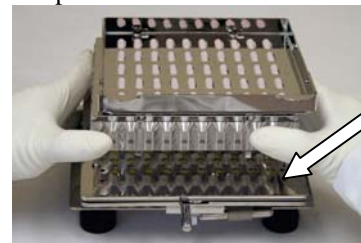

6. To fill remaining rows, repeat steps 1-4 except this time align the Orienter locating feet in the holes nearest the operator.



Use the Tablet Filler Attachment with the Profill 100 or Profill OE visit www.capsuleworld.com For more information

DESIGNED & MANUFACTURED BY

TORPAC®

www.torpac.com info@torpac.com
T: 1-973-244-1125 F: 1-973-244-1365
333 Route 46, Fairfield, NJ 07004,