







**The Quickest Way to do Tablet Over-Encapsulation.**


-  Simple to use.
-  Works with the ProFill 100 & ProFill OE Capsule Filling Systems.
-  4 Easy-to-change tablet sheets allow you to fill a variety of round tablets, from 6.1 to 9.0 mm in OD and 2.0 to 5.0 mm in height.
-  Custom tablet sheets are available for tablets less than 9.0 mm OD but outside the above size range

*The ProFill Capsule Filling System not included.*

**How to Over-Encapsulate**



**1.** Use Tablet Sheet selector on page 2 to determine the correct tablet sheet to install in the TFA.

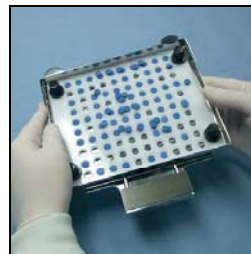

**2.** Pour 100+ tablets on to the TFA. Shake back and forth moving them into slots.



**3.** Lift gate and pour off excess tablets.



**4.** Place the TFA on ProFill System after outer capsules have been separated.



**5.** Press the release mechanism to drop the tablets into the capsules. Remove the TFA.

*Tip: You may wish to fill the capsules with a suitable excipient to avoid rattling.*



**6.** Replace the ProFill System Caps Tray and lock the capsules as per normal operation.



**Use the Tablet Filler Attachment with the Profill 100 or Profill OE visit [www.capsuleworld.com](http://www.capsuleworld.com) For more information**

**DESIGNED & MANUFACTURED BY**



www.torpac.com info@torpac.com  
T: 1-973-244-1125 F: 1-973-244-1365  
333 Route 46, Fairfield, NJ 07004, USA

## Tablet Sheet Selector

Recommended Tablet OD Range	Tablet Height Range		Tablet Sheet Label
	Recommended	Double Tableting may occur*	
6.1 to 7.0	2.5 to 3.4	2 to 2.4	<b>A</b>
7.1 to 8.0	2.5 to 3.4	2 to 2.4	<b>B</b>
8.1 to 9.0	2.5 to 3.4	2 to 2.4	<b>C</b>
7.6 to 9.0	3.5 to 5.0	2 to 3.4	<b>D</b>


For tablets less than 6.1 mm in OD or more than 2 mm thick, custom plates are available.  
\*In this height range you may get 2-3 tablets in a tablet sheet hole. This is easily noticeable and the extra tablets can be removed by hand.



Tablet OD Range

Tablet Height Range


## Changing Your TFA Tablet Sheet

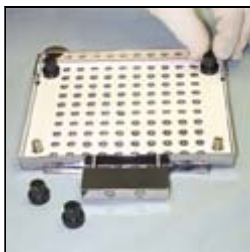


1. Remove the four thumbscrews from the TFA Base, and remove the white Tablet Sheet.



2. Select white Tablet Sheet using the Tablet Sheet selector above.

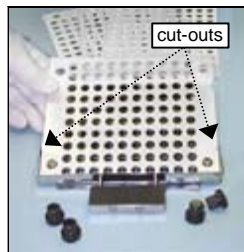

3. Install Tablet Sheet on the TFA base. Align the four corner holes with the screw posts on the TFA base and position the two cut outs to the front.

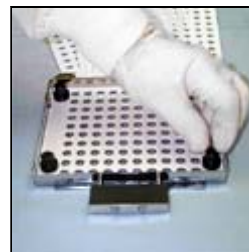

4. Replace the four thumbscrews removed in step 1. You are now ready to use your TFA for a new tablet size.

