



#9 / 9el Multi Capsule Filler

Operating Instructions

Ideal for filling multiple capsules at one time when fill weights can be within +/- 7 to 10% of average (e.g. pharmacokinetic studies)

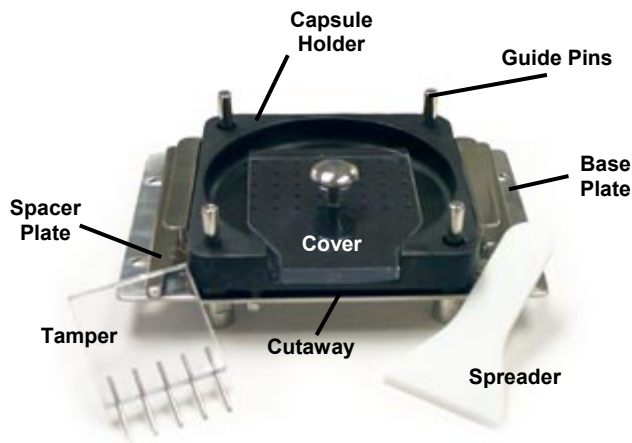


1. Pull gloves tight around fingertips for easier capsule handling.

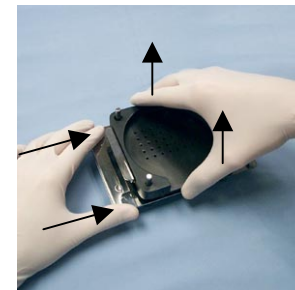


2. Lift the Capsule Holder and push the Spacer Plates in and under the Holder.




3. Separate capsules one at a time and place caps in a shallow container such as a Petri dish.

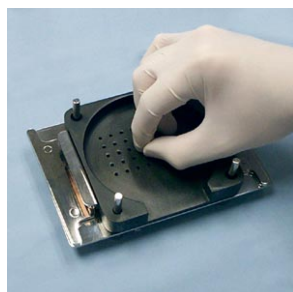


4. Place the capsule body (longer, bottom half) into the Capsule Holder with the body opening facing up. Do this for all 40 capsules.

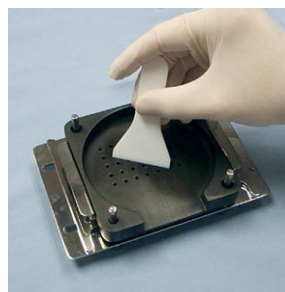


5. When all the holes have capsule bodies in them, pour out a pre-measured quantity of powder. Spread the powder with the Spreader.

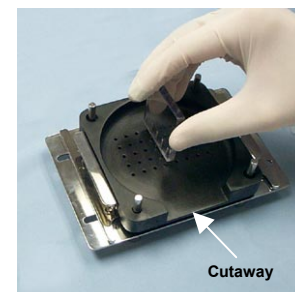

6. Use the Tamper to pack powder into capsules. Spread remaining powder as in step 5 for proper fill weight. Push the excess powder out through the cutaway into a weighing boat or other container.

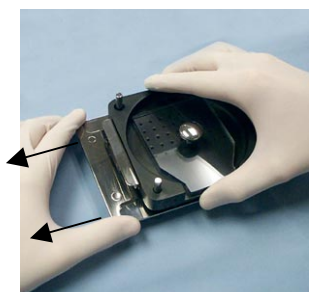


7. Place the Cover on top of the Holder. This ensures capsules stay fully seated. Now, hold the Capsule Holder and slowly pull out both Spacer Plates. Next, gently let down the Holder.

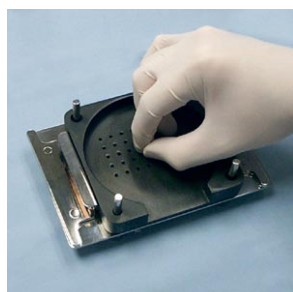

8. Remove the cover. Place the caps back onto the capsule bodies, pressing to fully lock each capsule.



9. Use the included Cover, pressing gently, to ensure the capsules are fully locked.

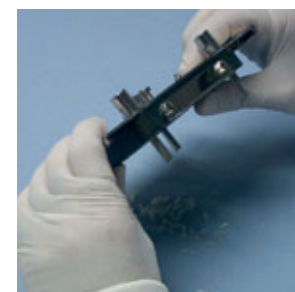

10. After all capsules are joined; turn the entire filler upside-down to remove all the filled capsules.

TORPAC[®]

www.torpac.com info@torpac.com
T: 1-973-244-1125 F: 1-973-244-1365
333 Route 46, Fairfield, NJ 07004, USA