Vet Clinic Filller

Sizes: Su07, 07, 10 (60-holes) up to 720 capsules/hour*

Size: 11 (70-holes)
up to 840 capsules/hour*

Sizes: 12el, 12, 13 (120-holes) up to 840 capsules/hour*

Sizes: 32150, 35110 (50-holes) up to 500 capsules/hour*





Watch video at www.torpac.com/vcf

For manufacturers a low-cost method to fill capsules for test-markets. For veterinary clinics a quick and easy way to fill capsules.

- Simple to operate, see video at www.torpac.com/vcf.
- Fills multiple capsules in one cycle. Manual orientation, separation & locking of capsules is required.
- Fill weights for individual capsules are within ±5% of average.
- Filler base includes base frame and 2 powder spreaders.
- Filler Change Parts include 2 filler plates and upper & lower spacers for size group.
- Easy to clean.
- Made from acrylic, nylon and lnox (stainless steel 316).
- Optional Tamper available as special order.



ProFil

Entry-level for R&D trials & clinics. Individually orient capsules. Use Filler to Separate, Fill, Tamp & Lock Capsules.

Sizes: Su7/7/10, 11, 12el/12/13 Output: 400+ Capsules/hour*



Capsule Veterinary System 2G

A bench-top system for regular production use. Approximately 3-5 million a year on 1 shift.

Sizes: Su7, 7, 10, 11, 12el, 12, 13 Output: 1000 - 2400 Capsules/hour*



AFV - Fully Automatic

Under Development.
Approximately 4-12 million capsule a year on 1 shift.

Sizes: Su7, 7, 10, 11, 12el, 12, 13 Output: 2500 - 7500 Capsules/hour*




ProMixer[™]10L

ProMixer Tumble Blender for mixing up to 10L (10 US Quart) of powder using V-Shell or Drum Shells. Typical mixing time 5-15 minutes based on % of air volume used and powders. Mixer & Shells made from Inox (Stainless Steel 316).

*Output based on powder, operator speed & number of taps/tamps.

Custom Dosing Applicator

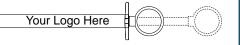

Designed & Manufactured by:

 $RPAC^{*}$

US - T: +1-973-244-1125 333 US 46, Fairfield, NJ 07004, USA

www.torpac.com info37@torpac.com

EU - T: +31-6-30-93-65-27 Jan Campertstraat 19, 6416 SG Heerlen, The Netherlands.

Available laser marked in black or embossed with your logo and product name. For use with sizes: 12el, 12, 13



Vet Clinic Filller

Size Su07-13 Capsules

These sizes use a common Filler Base. Order Change Parts for capsules sizes you want to fill.


Item Description		Capsule Size	Item #	US\$
Filler Base Kit Su07-13 (requires Change Parts) Base Kit includes: 1 Bottom Plate, 4 rubber feet, 4 Locating Pins, 4 Nuts, 2 Spacer Pins, 2 Upper Spacers & 2 Powder Spreaders (not shown in photo). Base Kit Su07-13 requires Change Part Set, see below.		Base Kit Su07 - 13	XV-VCF FB	\$220
Change Parts Set		Size Group		
Change Parts Set includes: 1 Top plate & 2 Lower Spacers common for an entire Size Group excepting 12el/12/13. Change Parts for sizes 12/12el/13 include a Top Plate and Lower spacers for 12el & 12. Size 13 requires unique Upper & Lower Spacers ordered separately for \$80. (XV-VCF FS 13) Required for use with Filler Base Kit Su07-13.		Su07, 7, 10	XV-VCF FCP A	\$450
		11	XV-VCF FCP B	\$375
		12el, 12, 13ª	XV-VCF FCP C	\$450
Tampers (optional) Optional Tampers to compress powder for increased fill weights & faster filling.		Su07, 07, 10	XV-VCF TMP A	\$190
		11	XV-VCF TMP B	\$200
		12el, 12, 13	XV-VCF TMP C	\$205

Size 32150 & 35110 Capsules


These sizes use a common Filler Base and are Matched with Charge Parts for a Size Group.

Item Description		Capsule Size	Item #	US\$
Filling Machine (Complete) The only filling system for Torpac King Size veterinary capsules. Includes 1 full system for either size 32150 or 35110.	and the same	32150	XV-VCF 32150	\$605
	- aminum	35110	XV-VCF 35110	\$605

Operating Overview

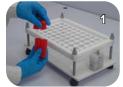

Separate Capsules

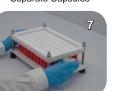

Rotate Upper Spacer



Insert Capsule Bodies



Replace Capsule Caps



Pour Pre-measured Powder



Capsules are Filled & Locked



Spread Powder



Tap or Tamp Powder



Spread remaining Powder

See Users Manual or Video for more Information at www.torpac.com/vcf