



Available in Sizes Su07/07/10, 11, 12e/12 in October 2012.

For manufacturers a low-cost alternative to an auger filler.  
For veterinary clinics a quick and easy way to fill capsules.

- Output per hour: Up to 500 capsule approx.
- Capacity: 50 capsules/cycle
- Simple to operate (see reverse)
- Fills Capsules. Manual orientation, separation & locking.
- Fill weights within  $\pm 5\%$  of average
- Includes everything you need to begin filling capsules
- Easy to clean.

| Veterinary Clinic Filler |              |       |
|--------------------------|--------------|-------|
| Capsule Size             | Item Code    | US\$  |
| 32150                    | XV-VCF 32150 | \$495 |
| 35110                    | XV-VCF 35110 | \$495 |

| Filling Veterinary Size Capsules  |   | Powder Mixing   | Filling Size 000-5 Capsules  |   |
|---|---|---|--|---|
|  |  |  |  |  |
| Hand-held Filler<br>ProFill Vet   | Bench-top filler<br>CVS   | 0.4 ltr / 0.8 Pint<br>Mixer / V-Blender   | Hand-held Filler<br>ProFill 100  | Bench-top Filler<br>ProFiller 3700  |

**Quick Start Guide**

Photos show filler for 30 capsules. Current model is filler for 50 capsules.



**1.** Separate the caps and bodies of capsules. Keep the caps in a bowl.



**2.** Place the bodies (longer half) in the filler, open end of the body up.



**3.** Pour pre-measured powder on to the Filler. When pouring, distribute powder evenly (*Tip A*).



**4.** Spread the powder evenly with the Powder Spreader.




**5.** While holding the top sheet and body sheet together lift the filler from one side by 1 cm (1/2 inch) and then drop it back on the table. Repeat for each of the four sides (one tapping cycle). See *Tip B*.




**6.** Spread any extra powder. Repeat tapping till all the powder is filled in the bodies. Typically, after 3-4 tapping cycles no more powder will enter bodies.

To fill additional powder into the capsule bodies, tap the filler (see next step).

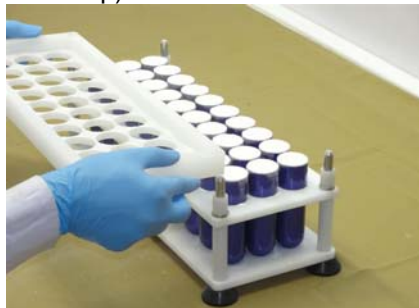

**7.** Remove the Top body sheet slowly.

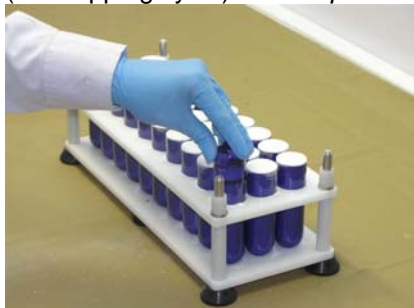

**8.** Place the caps back on the capsule bodies. Press down to ensure the cap locks on to the body (*Tip C*).




**9.** Remove capsules from the filler by turning it over. With practice you can fill up to 500 capsules/hour.

**Tip A:** If it appears that powder will flow over the top body sheet stop and go to step 4. When there is space on the Top Body Sheet add remaining powder.

**Tip B:** Video demonstrating tapping is at [www.torpac.com/vcf](http://www.torpac.com/vcf).

**Tip C:** When locking the first resistance you feel is the lock ring on the cap stops at the top of the body. Push down firmly until the lock ring on the cap engages with the matching lock ring on the body.

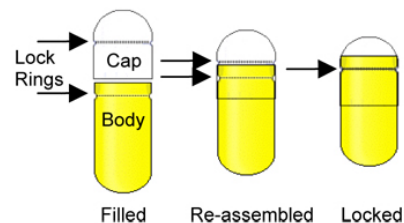

**Cleaning:** Made from acrylic. Hand-wash with soap & water. Do not clean in a dishwasher or use solvents.